

# Sciencescope *TrueColor* CaptureX

## Installation Guide

### **Windows XP**

Make sure the camera **isn't** connected to the computer

**Run** setup file and follow prompts \*

**Click** finish

**Plug in** the camera to a USB 2.0 port

Windows will prompt you with a hardware installation wizard

Let Windows search for the drivers automatically \*\*

**Click** Finish

**Run** CaptureX from either the desktop or start menu \*\*\*

### **Windows Vista**

Make sure the camera isn't connected to the computer

**Right click** on the setup file and run as administrator

Follow prompts

**Click** finish

**Plug in** the camera to a USB 2.0 port

Windows will prompt you with a hardware installation wizard

Let Windows search for the drivers automatically \*\*

**Click** Finish

**Run** CaptureX from either the desktop or start menu \*\*\*

\*Windows XP Service Pack 3 requires workaround, please see WinXP SP3 workaround guide

\*\*If Windows was unable to locate the drivers, please see troubleshooting guide for manual installation

\*\*\*If you encounter a black screen, please see troubleshooting guide on how to reset video settings